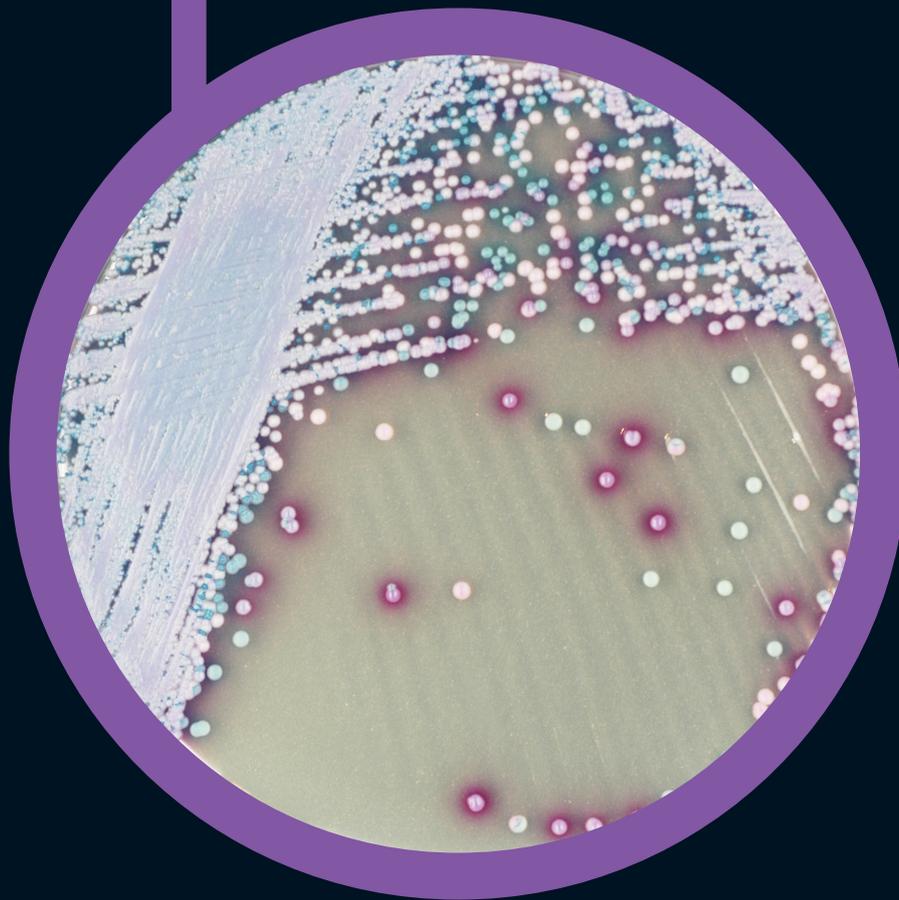
CHROMagar™ Malassezia

For detection of *Malassezia* spp.

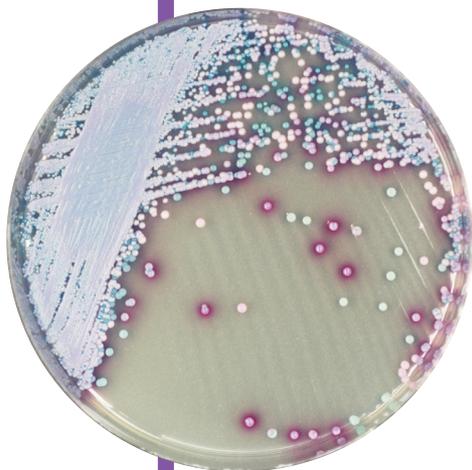
Plate Reading

For detection of:

- *Malassezia furfur*
→ large, pale pink and wrinkled
- Other *Malassezia spp*
including *M. globosa* & *M. restricta*
→ mostly pink to purple

For detection of *Malassezia spp.*

Background

Malassezia is a fungi naturally found on the skin of many animals, including humans. On normal healthy skin it does not cause infections, but when the environment of the skin is altered, *Malassezia* species are able to cause several cutaneous diseases as severe dermatitis or otitis (inflammation of the skin or ears respectively)

Medium Performance

1

EASY AND COLOURFUL DETECTION

Since members of the genus *Malassezia* share similar morphological and biochemical characteristics, the use of traditional culture media for differentiating them based on phenotypic features is not suitable. CHROMagar Malassezia was developed with the goal of facilitating not only their detection, but also to improve an algorithm for the differentiation of the most common species (see: Revised Culture-Based System for Identification of *Malassezia* Species, by Takamasa et al. JCM No-2007)

Typical Appearance of the colonies reported in the study:

<i>M. pachydermatis</i> CBS 1879 → large, pale pink & smooth
<i>M. restricta</i> CBS 7877 → small, pink & smooth
<i>M. dermatitis</i> JCM11348 and JCM11470 → small, pale pink & smooth
<i>M. slooffiae</i> CBS 7956 → large, pale pink & smooth
<i>M. obtusa</i> CBS 7876 → medium, pink & rough
<i>M. globosa</i> CBS 7966 → small, purple & smooth
<i>M. sympodialis</i> CBS 7222 → large, pale pink & smooth
<i>M. furfur</i> CBS 1878 → large, pale pink & wrinkled

Extracted from: «Revised Culture-Based System for Identification of *Malassezia* Species», by Takamasa et al. JCM No-2007)

Medium Description


Powder Base	
Total	56.3 g/L
Agar	15.0
Peptones and extracts	38.0
Chromogenic mix	2.8
Chloramphenicol	0.5
Storage at 15/30°C - pH: 5.8 +/-0.2	
Shelf Life	3 years

Mandatory Supplements to the Powder Base (Not provided by CHROMagar):

Glycerol	2 g
Tween 40	10 g

Usual Samples	Clinical and veterinary samples (skin, ear canals etc.)
Procedure	Direct Streaking. Incubation 72h at 30-37°C Aerobic conditions.

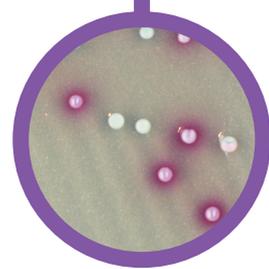Scientific Publications on this product: available on www.CHROMagar.com
Please read carefully the instructions for use (IFU document) available on www.CHROMagar.com

Quality Control Strains

<i>M. furfur</i> ATCC® 14521	pale pink
<i>C. albicans</i> ATCC® 10231	pale green
<i>C. tropicalis</i> ATCC® 1369	steel blue + mauve halo
<i>C. glabrata</i> ATCC® 2001	mauve
<i>C. krusei</i> ATCC® 14243	pink
<i>E. coli</i> ATCC® 25922	inhibited

From Takamasa et al. study:

<i>M. pachydermatis</i> CBS 1879 → large, pale pink & smooth
<i>M. dermatitis</i> JCM11348 and JCM11470 → small, pale pink & smooth
<i>M. globosa</i> CBS 7966 → small, purple & smooth

ATCC® is a registered trademark of the American Type Culture Collection


Order References

Please use these product references when contacting your local distributor:


5000 ml pack	MZ282
25 L pack	MZ283-25

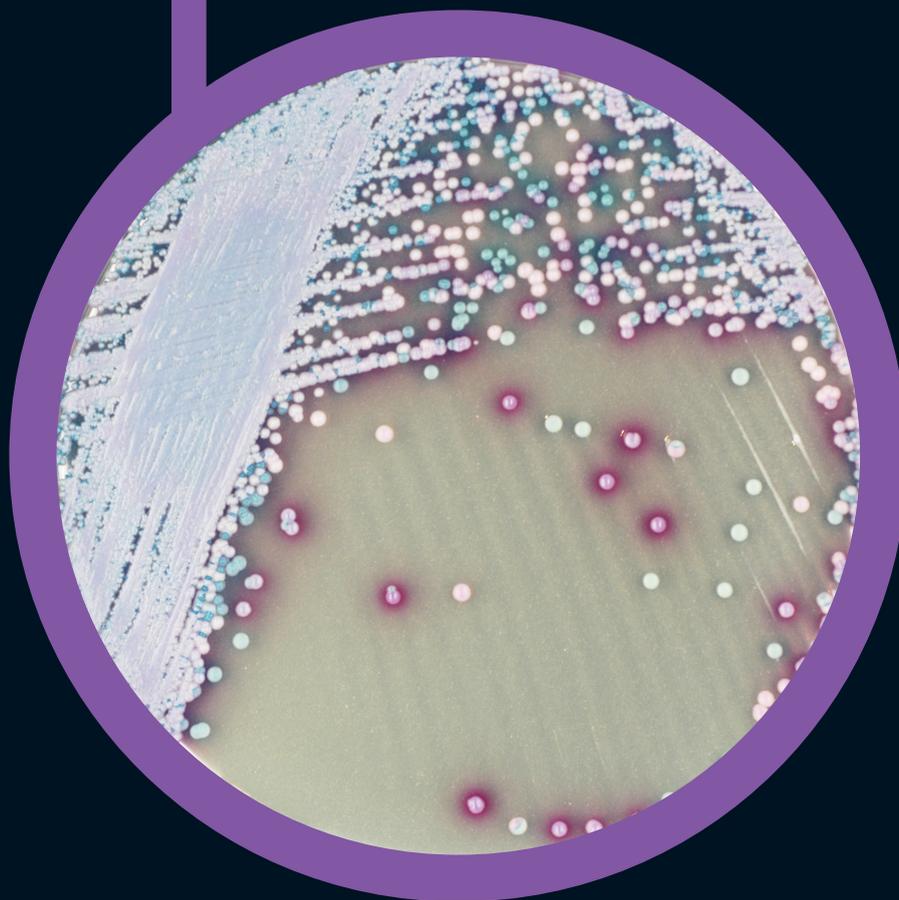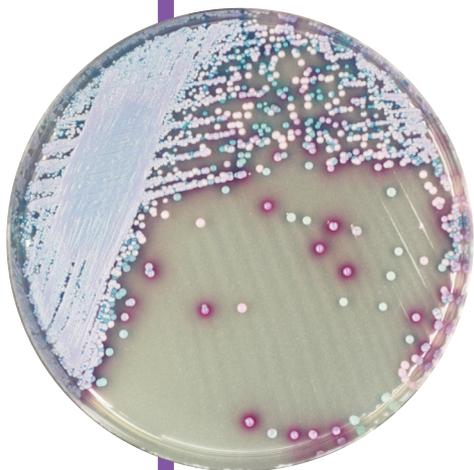Manufacturer: CHROMagar
4 place du 18 juin 1940 75006 Paris - France
Email: CHROMagar@CHROMagar.com
Website: www.CHROMagar.com
Find your nearest distributor on
www.CHROMagar.com/contact

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